QUICK  
START  
GUIDE

**Moticam X<sup>3</sup>**

EN

### Quick Start Guide

Thank you for purchasing this Moticam. This camera is a delicate instrument and should be handled with care. When not in use, please store the camera in a safe place with the dust cap attached. Do not let the camera get in contact with water or expose it to extreme temperatures (below 0 or above 50 degrees Celsius). In order to receive the highest level of service, please retain your receipt as proof of purchase and **register your camera online at [www.motic.com](http://www.motic.com) today.**

Should you have any questions please do not hesitate to **contact our customer service.**

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**EN This device conforms to US FCC and european EMC regulations:** operation is subject to the following two conditions:

1. This device may not cause harmful interference, and
2. This device must accept any interference received, including interference that may cause undesired operation.

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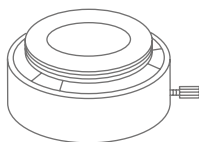
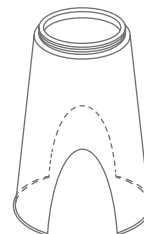
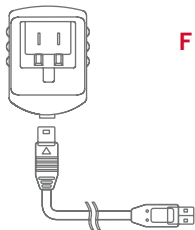
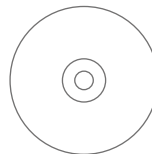
### EN Using the camera on an eyepiece

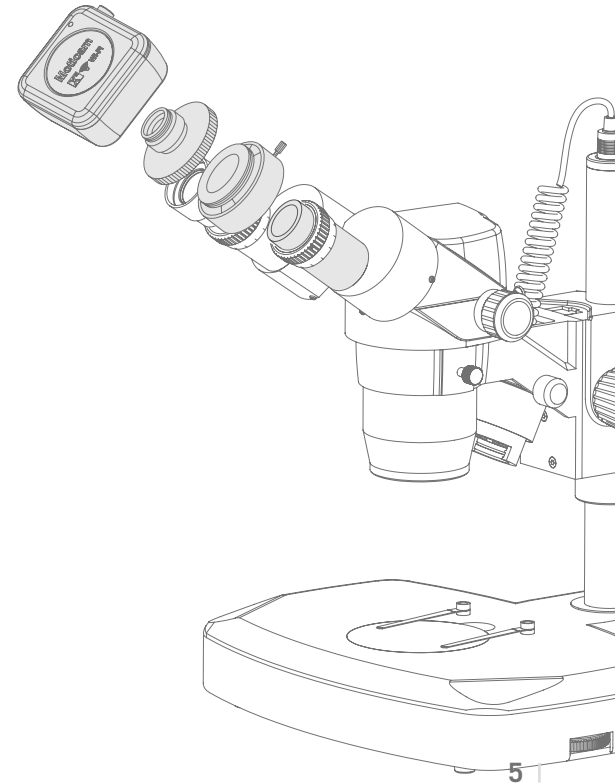
1. Remove the Dust Cap (B) from the camera (A) and attach the focusable lens (D).
2. Choose the right size eyepiece adapter (E) and attach this to the focusable lens.
3. Attach this assembly to your microscope's eyepiece by tightening the centering screws on the eyepiece adapter.
4. Connect the USB power supply to the camera. When the light on the camera is on, it means the camera is working and transmitting images and you can connect to it through WiFi.
5. Connect to the Moticam X<sup>3</sup> WiFi signal. Please see the instructions on page 11.

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### EN Using the camera as a Macro Camera

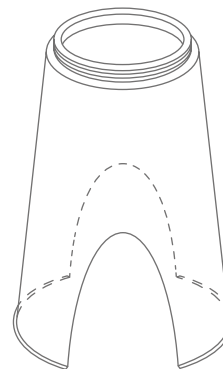
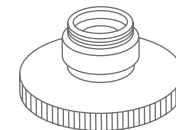
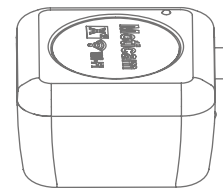
1. Remove the Dust Cap (B) from the camera (A) and attach the focusable lens (D).
2. Attach the Macro Tube (H) onto the lens
3. Place this assembly over the item you wish to view and ensure there is adequate lighting. If necessary, a desk lamp is a good solution.
4. Connect the USB power supply to the camera. When the light on the camera is on, it means the camera is working and transmitting images and you can connect to it through WiFi.
5. Connect to the Moticam X<sup>3</sup> WiFi signal. Please see the instructions on page 11.

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### EN Using the camera on a Trinocular Microscope

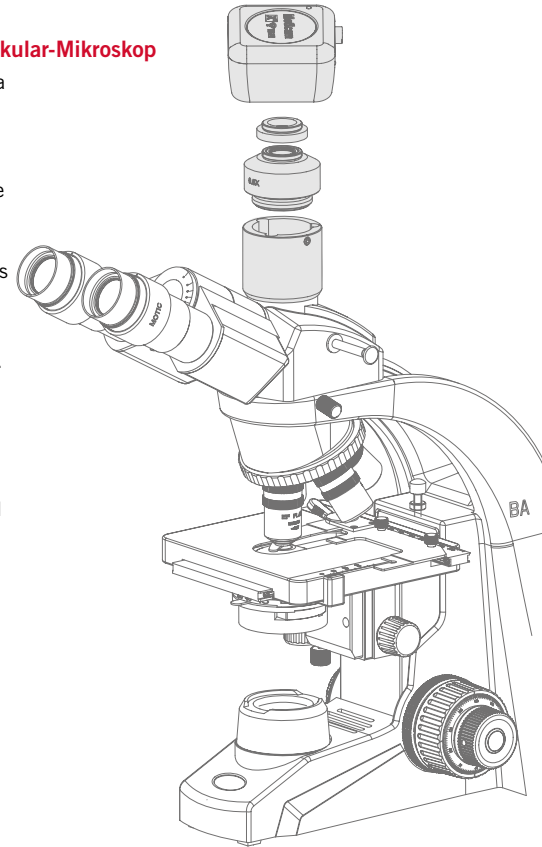
1. Remove the Dust Cap (B) from the camera (A) and attach the microscope's C-Mount adapter (\*Please note that the C-Mount adapter is specific to your microscope. For Motic microscopes, please consult the relevant Motic Accessory list and choose an appropriate adapter for your Moticam. Depending on the type of adapter and microscope, the C-Ring (C) may be necessary between the camera and the C-mount adapter).
2. Connect the USB power supply to the camera. When the light on the camera is on, it means the camera is working and transmitting images and you can connect to it through WiFi.
3. Connect to the Moticam X<sup>3</sup> WiFi signal. Please see the instructions on page 11.

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### **EN Connect to the Moticam X<sup>3</sup> WiFi signal**

1. Each Moticam X<sup>3</sup> transmits its own WiFi signal to which you can directly connect your WiFi Device.
2. The Moticam X's SSID signal name is printed on a label on the camera. To log into your camera, simply connect your WiFi device to that signal. The password for the network by default is 12345678.

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## **EN Using a Wifi device that operates with Android or iOS**

\* For best performance, we recommend to use WiFi devices, such as tablets, with a screen not smaller than 7 inches.

For iOS devices we suggest using iOS 5 or higher while for Android Devices we suggest using Version 4.0 or higher.

1. To operate and use the camera on your tablet, please download our free App called MotiConnect. For iPad you can find this on AppStore, for Android devices you can find this on Google Play.
2. You can also use certain web browsers such as Chrome, Firefox or Safari to view, measure and capture without the need of the App.  
To view the camera stream on your browser, please type in the IP number also shown on the label on the camera. For example 192.168.1.151.  
On your browser, type 192.168.1.151:8080 and you will be connected straight to the camera.
3. The camera supports both MJPG and H.264 output. By default the camera is set to H.264 in order to provide fast FPS transmission to the MotiConnect App.  
If you would like to view the live image on both MotiConnect and a Web browser, please change the output mode to MJPG using the settings on the MotiConnect App.

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### **EN Use the Moticam X<sup>3</sup> on your Wifi enabled computer**

The Moticam X<sup>3</sup> can also be used with Motic Images Plus software for computers. If you have a WiFi enabled laptop, you can go to [www.motic.com](http://www.motic.com) and download Motic Images Plus 3.0ML software by registering as a user with the DM serial number you will find on the box. To view the camera in our product's MI Devices interface, simply log your computer onto the Moticam X's WiFi signal, by selecting this Wifi connection and entering the password. Then open the MI Devices interface by clicking on the camera icon within Motic Images Plus and select Moticam X<sup>3</sup> from the drop down menu. Then click on Open.

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