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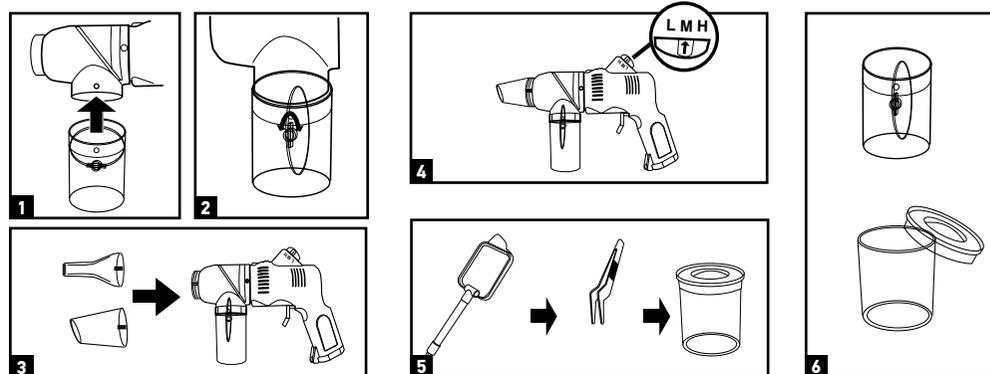
Levenhuk, Inc. 924-D East 124th Ave. Tampa, FL 33612 USA  
Levenhuk®, LabZZ® are registered trademarks of Levenhuk, Inc.  
© 2018 Levenhuk, Inc. All rights reserved.  
[www.levenhuk.com](http://www.levenhuk.com)  
20180726

## EN Levenhuk LabZZ BC10 Bug Collector

**May be used by kids over 3 years old. Observe insects through the container only, do not touch them – some insects can bite or sting you. Make sure to open the container and release the insect when your observation is finished.**

Levenhuk LabZZ BC10 Bug Collector

Levenhuk LabZZ BC10 Bug Collector is an excellent kit for fun and safe collecting and studying spiders, flies, grasshoppers, mosquitos, caterpillars, butterflies, and many other insects.

**Using:**

- Install the container as shown (1). Rotate the cover of the container to open it (2). Choose and install one of the nozzles (3).
- Place the nozzle next to the bug and press the vacuum button. The Bug Collector is equipped with three speed levels: „L” is for low speed, “M” is for medium speed, and “H” is for high speed. Select the speed depending on the size of the insect (4).
- You can also use a scoop-net or tweezers to catch insects (5).
- Open the cover of the container or a glass to release the insect (6).

**Batteries installation and replacement:** use a screwdriver to open the battery compartment cover, install 4 AA batteries and close the cover. A screwdriver and batteries are not included in the kit.

**⚠ Always replace the whole set of batteries at one time; taking care not to mix old and new ones, or batteries of different types. Remove used batteries promptly. Never heat batteries in order to revive them. **Seek medical advice immediately if a small part or a battery is swallowed.****

Warranty: 6 months. For further details, please visit our web site: www.levenhuk.com/warranty
The manufacturer reserves the right to make changes to the product range and specifications without prior notice.

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